USB Interface Kit
P/N 71208792
Instruction Manual

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1. INTRODUCTION

This USB Interface Kit P/N 71208792 provides a solution to connect an Adventurer Pro balance to a computer via a Universal Serial Bus (USB) port. This interface kit uses a USB-serial conversion class that allows the use of a USB port like an RS232 serial port, also known as a Virtual COM port. Please read this manual completely before installation and operation.

READ ALL SAFETY WARNINGS BEFORE INSTALLING, MAKING CONNECTIONS, OR SERVICING THE BALANCE. FAILURE TO COMPLY WITH THESE WARNINGS COULD RESULT IN PERSONAL INJURY AND/OR PROPERTY DAMAGE. RETAIN ALL INSTRUCTIONS FOR FUTURE REFERENCE.

1.1 Contents

- USB Printed Circuit Board (PCB)
- PCB Cable
- USB Cable
- Protective Cone Cover
- Screws (2) to secure the USB PCB
- Program CD

1.2 Tools and PC System Requirements

- Philips screwdriver or Torx driver T10 & T20 (depending on model)
- PC USB port
- Windows 98[®], Windows 98SE[®], Windows ME[®], Windows 2000[®] or Windows XP[®]

2. INSTALLATION OF HARDWARE

Install the hardware first before the USB driver software contained in the kit program CD.

2.1 Safety Precautions

- Disconnect the balance from the mains power supply and remove any installed batteries before beginning installation.
- Internal components, such as the load cell and printed circuit boards (PCB), are delicate and need to be handled with care.
- Do not subject the internal components to static electricity discharge.
- The PCB's should be handled by grasping the edges only and never placing fingers on any of the runs or traces.

2.2 Hardware Installation

The Adventurer Pro series comes in different configurations (see Figures 1A-B-C). Disassemble/assemble the balance according to the configuration that matches your model.

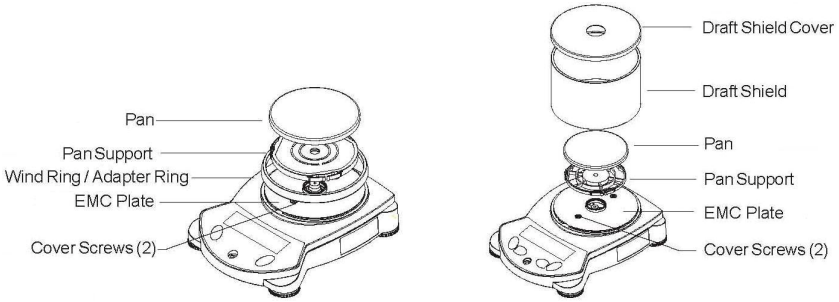
Figure 1A.

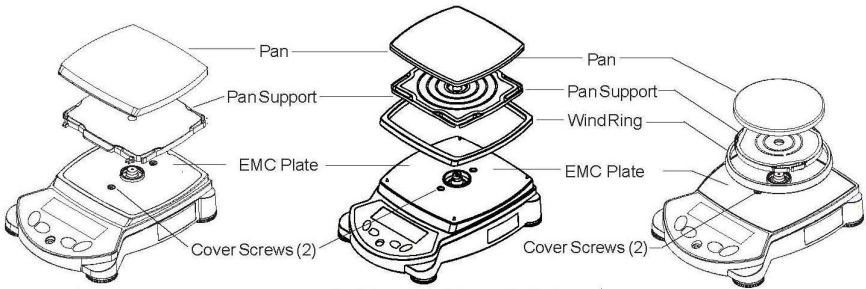
Figure 1B.

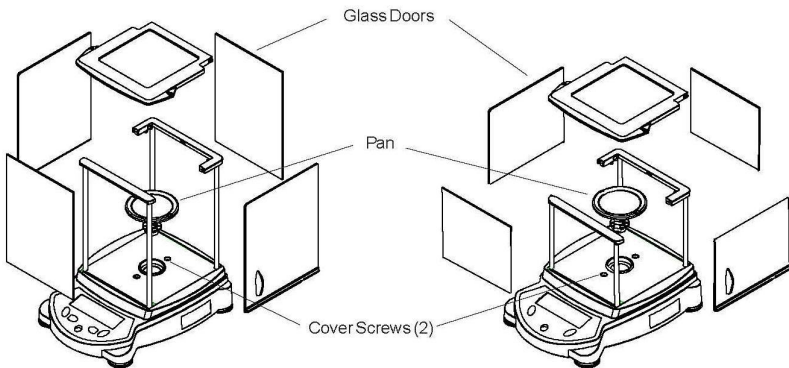
Figure 1C.

1. With power removed from the balance, remove any draft shield or draft shield glass panels, the pan, pan support and any wind ring or adapter ring (sizes and shapes vary) to fully expose the EMC plate on the top of the balance.
2. Remove the 2 screws and the EMC plate.
NOTE: On some models with angular wind rings, be careful not to lose the 4 corner plastic locating pins where the EMC plate slots over. These are marked “F” (front) or “B” (back) with the markings facing inwards (note their positions in case dislodged and needing to be re-installed).
3. Disassemble the top from the bottom housing (depending on model):
 - a) Small frame models: Remove the 4 feet, then remove the 4 screws inside the openings uncovered by the removal of each foot.
NOTE: While removing the feet, avoid resting the balance on its top side (over the load cell cone assembly). If needing to do so, place the protective plastic cone cover from the kit over the load cell cone assembly to minimize risk of damage to the load cell (See Figure 2A).
 - b) Large frame models: Remove the 2 screws (and any grounding clip) on the top of the balance that were exposed by the removal of the EMC plate.
NOTE: If the balance was trade approved, any sealing sticker covering the housing screws will need to be removed. In effect, the balance will have to be re-approved and re-sealed.
4. Carefully lift the top cover from the bottom housing, mindful of the cable connected to the main PCB assembly.
5. Unplug the cable from the main PCB assembly, then set aside the top cover.
6. Remove the 2 screws securing the COM2 cover plate towards the back and right of the bottom housing (See Figure 2B). Discard the cover plate and 2 screws.
7. Mount the USB PCB over the corresponding stands, then secure with the 2 screws provided (See Figure 2C).
8. Connect the PCB cable to: a) the USB PCB, and b) to connector CN6 on the main PC board on the front of the balance (see Figure 2D). Route the PCB cable towards the left side of the bottom housing.
9. Re-assemble the balance in the reverse order of the disassembly process, being careful not to pinch any cables within or between the top and bottom housing.

Cone Cover (for small-frame models)

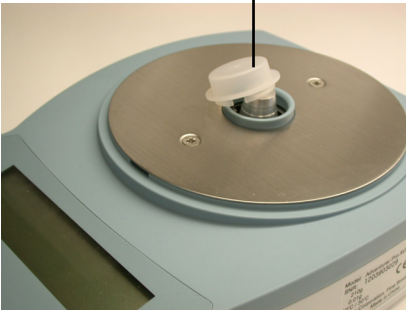
Figure 2A.

COM2 Cover Plate

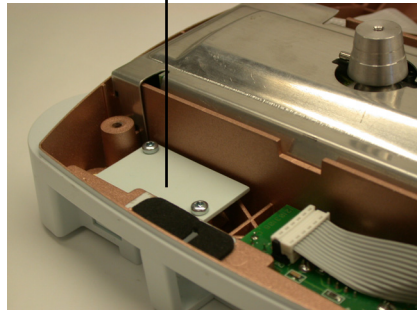
Figure 2B.

USB PCB

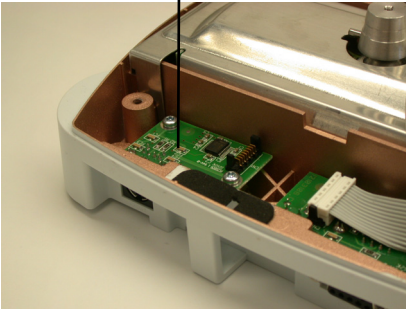
Figure 2C.

PCB Cable


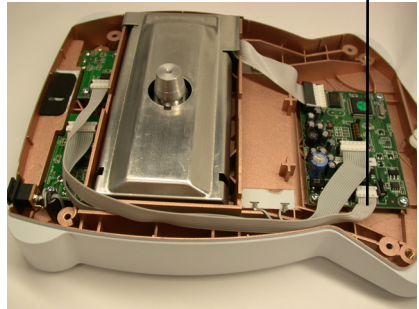
Figure 2D.

3. INSTALLATION OF SOFTWARE

After installing the hardware, install the USB driver software contained in the kit CD.


3.1 Software Installation

1. Insert the supplied CD into the CD-ROM drive of the computer.
2. Connect the balance to the computer's USB port through the USB cable (supplied).
3. "Found New Hardware Wizard" window will come up. The following screen shots are examples of the installation using Windows 2000®. Other Microsoft Windows® operating systems may show a similar window with a slightly different look. "Hardware Wizard" in all versions of Windows® will guide you through the installation of the necessary 2 drivers located on the CD with steps described below.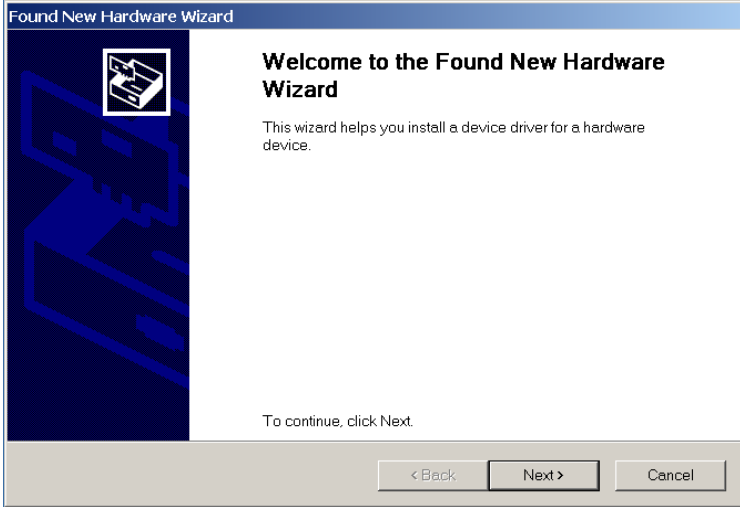


SCREEN 1

4. Click **Next** and screen 2 will be displayed. Click **Next** to move to screen 3 and check **CD-ROM Drive** before clicking **Next**.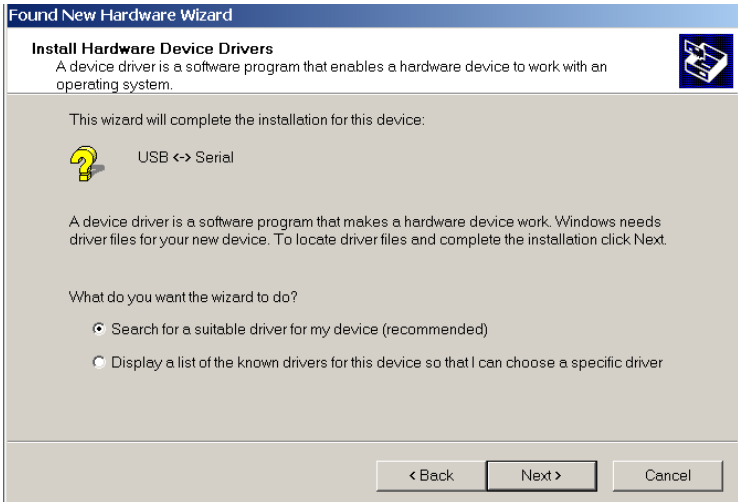


SCREEN 2

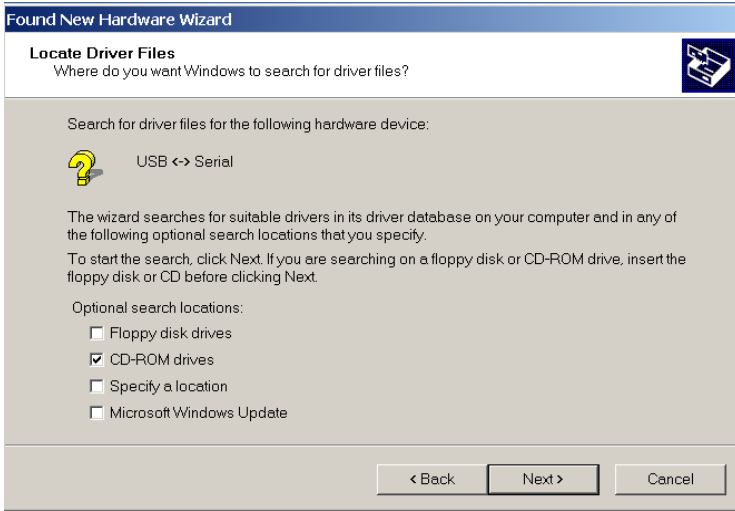SCREEN 3

- Two drivers will be installed in back-to-back sequence. Therefore, you'll repeat the steps in screen 1 to screen 4 to complete installation.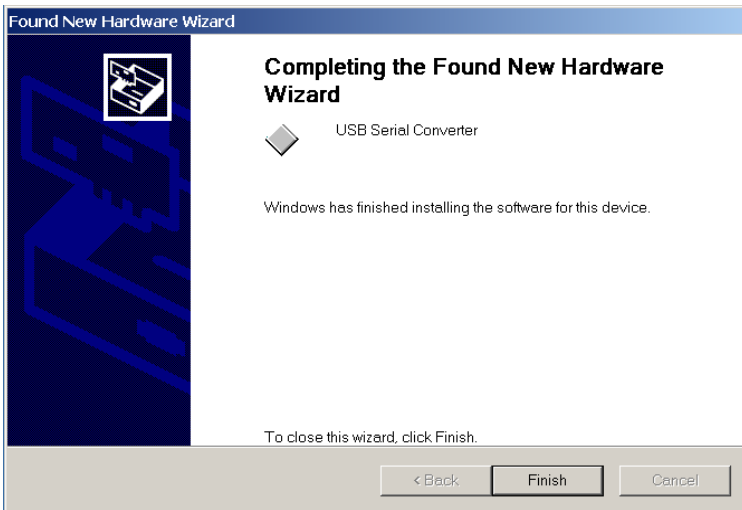


SCREEN 4


- After clicking the second **“Finish”** button, the USB driver software is installed and Windows® assigns a Virtual COM Port to the USB interface. Windows® typically assigns COM Ports in number sequence after previous assignments. For example, if COM3 and COM4 have been previously assigned, the Virtual COM Port for the new USB interface (which is also named USB Serial Port in Windows®) will be associated with COM5.

3.2 Finding USB Serial Port Information

Information on the USB Serial Port can be found by going to the control panel and clicking on **“SYSTEM”**. A System Properties window will come up. Click the **“Hardware”** tab and click the **“Device Manager”** button to bring up the Device Manager window.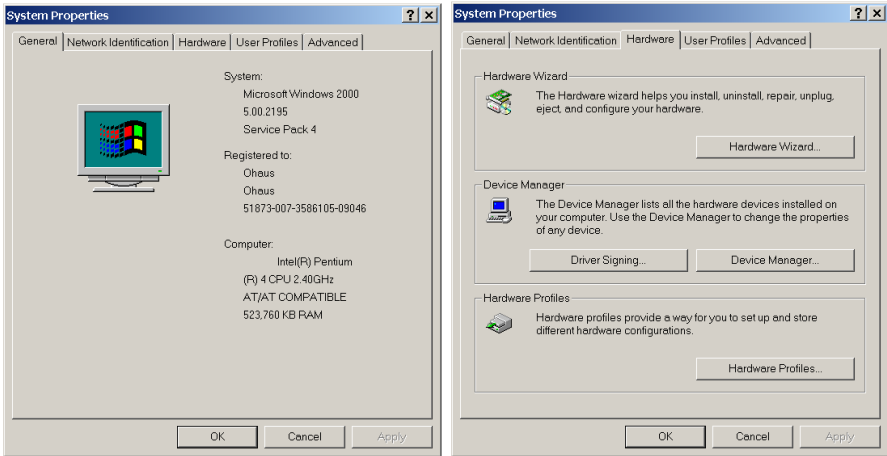

The USB Serial Port can be found under Ports (COM & LPT) in the Device Manger window. In the case of the picture shown below, the Device Manger window shows the COM port number of the USB Serial Port as COM3. The COM port number shown in this window is the COM port number the PC application will use to communicate with balance.

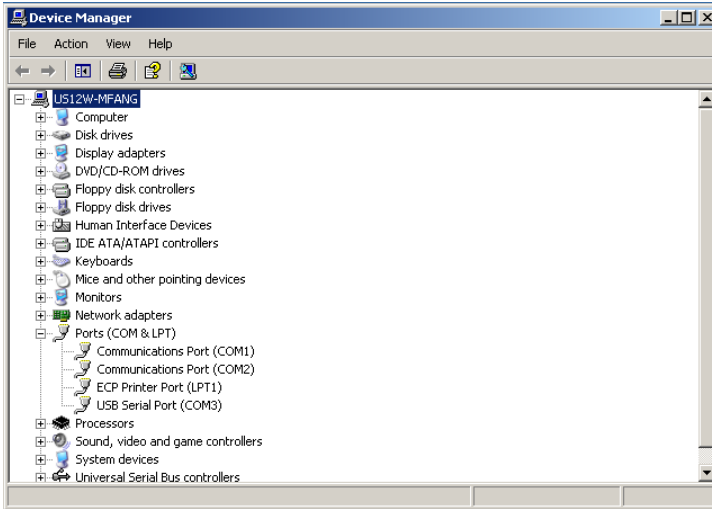Note: When using the USB interface with programs that limit the number of COM port designations (e.g. Ohaus MassTracker allows only COM1, 2, 3, & 4), it may be necessary to assign one of these port numbers to the new virtual port. To change the port number assigned to USB Serial Port, double click “**USB Serial Port**” in the Device Manager window. A “USB Serial Port (COMx) Properties” window will come up. Click the “**Port Settings**” tab and click on the “**Advanced**” button to bring up “Advanced Settings for COMx” and change COM port number.

4. BALANCE MENU SETUP

The Adventurer Pro software defines the USB Serial port as the RS232-2 port. Refer to the instruction manual supplied with the Adventurer Pro balance. Set up the RS232 communication and printing parameters as described in the RS232-2 and Print-2 submenus (enabled when this kit is installed).

5. MAINTENANCE

Service Information

For service assistance in the United States, call toll-free 1-800-526-0659 between 8:00 AM and 5:00 PM Eastern Standard Time. An Ohaus Product Service Specialist will be available to assist you. Outside the USA, please visit our website at www.ohaus.com to locate the Ohaus office nearest you.

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P/N 80100401 A © Ohaus Corporation 2010, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati.
Ohaus Corporation