



# 1 Create Account

Visit <https://www.traceablelive.com>  
Click on New Account.

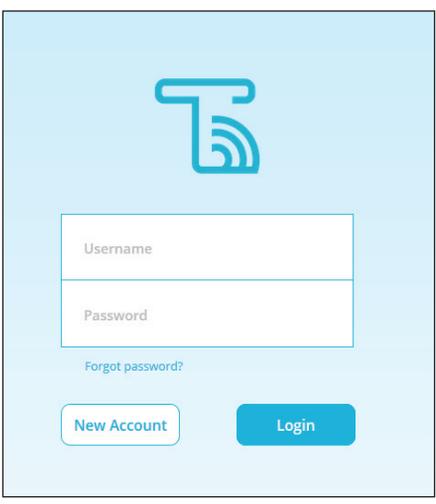


# 2 Enter Account/User Information

Fill in all required fields.  
**Note:** Password must be a minimum of 8 characters including a number and a special character (!@#%&^\*).

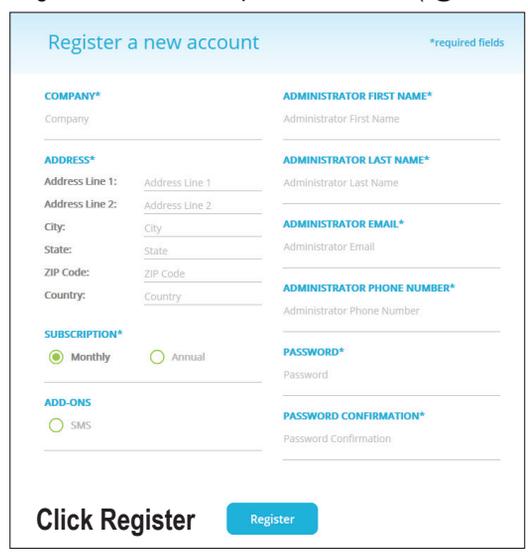

# 3 Enter Payment Info./ Confirm Charge

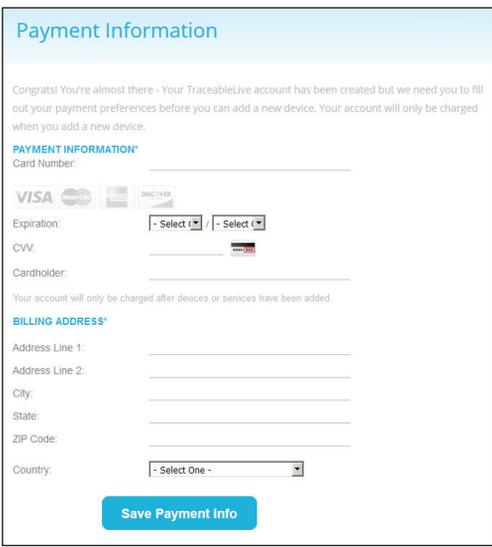

If this screen does not appear, go to log in screen, account settings and add payment information.  
**Note:** Payment must be set up prior to adding a device to your service.  
**Click Save Payment Info**

# 4 Begin WiFi Setup



On device Press/HOLD WiFi button for 3 secs. "RP" will appear on display.


Wait for a few seconds then press (no need to hold) WiFi button, "RP READY" will appear on display.




5 Commonwealth Ave  
Woburn, MA 01801  
Phone 781-665-1400  
Toll Free 1-800-517-8431

# Connect Unit to WiFi network via Mobile device

5

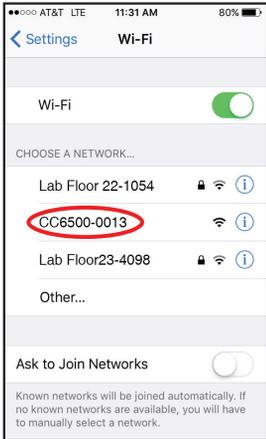

Using a WiFi enabled device (smart phone, tablet, etc.), find **CC6500-XXXX** in available networks. XXXX will be last 4 digits of unit's S/N# (on back of unit).

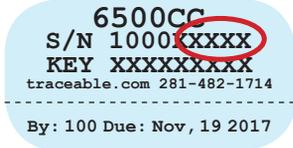

# Open Web Browser on your mobile device

6



Enter into web browser: **192.168.1.1**

Select desired local network to which the device will be connected, and enter password for network.

The most common network is **WPA2**.

# Press Add/Confirm device restart

7

Click Add



Unit will reboot, connect to WiFi, and begin posting data.

**Note:** Upon successful connection, you will receive an error message in your browser (Example: Safari cannot open the page because the network connection was lost.), and device reboot may take up to 1 minute, mobile will exit WiFi.

**Note:** The WiFi symbol will flash on device display until the device posts data successfully.

**Note:** ALLOW UP TO 30 MINS FOR DEVICE DATA TO UPDATE AND START TRANSMITTING DATA.

If and ERR message occurs on device, press acknowledge button (✓) and repeat steps 4-7.


Log-in to TraceableLIVE account to manage device settings, alarms, users, and payment options.

# Enter Device Information (located on back of unit)

8

Go to [www.traceablelive.com](http://www.traceablelive.com)

Click on Add New Device

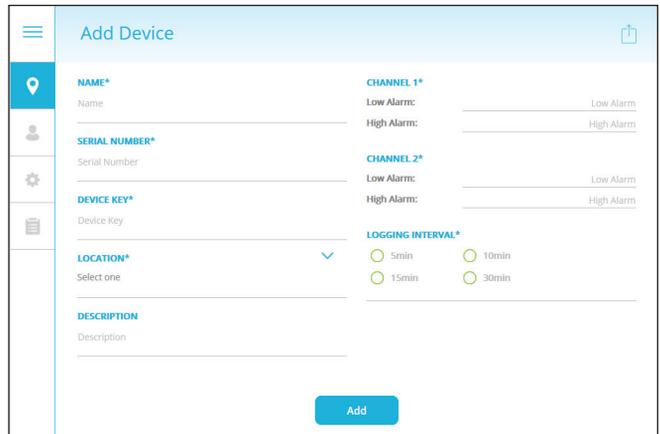

Complete all required fields.

Enter S/N & Device Key

Device and account setup is now complete.

